

Hotfix 4.17.1.3-1.1

Subject of Changes

Hotfix 4.17.1.3-1.1 fixes the following problem(s):

- When opening the reagent kit status window ("Functions" / "Reagent kit status"), it takes approx 1 minute till the window opens. The same occurs if the user presses one of the "Set full", "Empty", "Replace all containers" buttons.

This was realized by changing a stored procedure in the database.

Changes compared to latest patch 4.17.1.3-1:

- BusinessObjects.dll of patch 4.17.1.3 is used
- The job file "Hotfix 4.17.1.3-1.1.xml" was changed and will now copy ConnectionString.txt and pw-cdx from the Cedex 2 directory to the hotfix directory.

General Information

- It is possible to de-install the hotfix
- It is possible to run the installation of the hotfix several times
- No conflicts with other hotfixes
- No data base limitation

Limitations

Limited to data base version 4.16 and 4.17

Installation

1. Execute self-extracting file „Hotfix 4.17.1.3-1.1.exe“
2. Select Cedex 2 directory (default: „C:\Cedex2“; use browse button to select different directory)
3. After extraction open subdirectory „Hotfix 4.17.1.3-1.1“ in the Cedex 2 directory (use windows explorer)
4. If Cedex 2 directory is not „C:\Cedex2“ open file „install Hotfix 4.17.1.3-1.1.bat“ with editor and change Cedex 2 directory path.
5. If Cedex 2 directory is not „C:\Cedex2“ open file „install Hotfix 4.17.1.3-1.1.xml“ with editor and change Cedex 2 directory path.
6. Install hotfix by double clicking “install Hotfix 4.17.1.3-1.1.bat”
7. Follow instructions displayed in the window “Installation Cedex 2 Hotfix 4.17.1.3-1.1”

De-Installation

It is not possible to de-install the hotfix.

Validation

This hotfix does not influence the analysis process and has no impact on measured and calculated results!

Cedex 2 **HOTFIX**



Test

The installation of this hotfix was successfully tested with the following Cedex 2 software versions:

- Cedex 2.1.1 Build 4.16.1.2
- Cedex 2.1.1.Build 4.17.1.3

Support

For further questions regarding the hotfix and its installation please contact:
support@innovatis.com

~ 082007.pm/ohe ~