

Hotfix 4.17.1.3-3

Subject of Changes

Hotfix 4.17.1.3-3 fixes the following problem(s):

- Takes the size of the new container for Cedex 2 standard systems (60mL each for Trypan Blue, cleaning solution, and detergent) into account.

This was realized by changing container size and liquid level in the data base. The values of Trypan Blue, Cleaning Solution and detergent are 60 (size) and 0 (liquid level). Therefore, it is necessary to change the reagent kit afterwards.

General Information

- It is possible to de-install the hotfix
- It is possible to run the installation of the hotfix several times
- No conflicts with other hotfixes
- No data base limitation

Limitations

Limited to data base version 4.13 to 4.20

Installation

1. Exit Cedex 2 GUI and Server application
2. Execute self-extracting file „Hotfix 4.17.1.3-3.exe“
3. Select Cedex 2 directory (default: „C:\Cedex2“; use browse button to select different directory)
4. After extraction open subdirectory „Hotfix 4.17.1.3-3“ in the Cedex 2 directory (use windows explorer)
5. If Cedex 2 directory is not „C:\Cedex2“ open file „install Hotfix 4.17.1.3-3.bat“ with editor and change Cedex 2 directory path.
6. If Cedex 2 directory is not „C:\Cedex2“ open file „install Hotfix 4.17.1.3-3.xml“ with editor and change Cedex 2 directory path.
7. Install hotfix by double clicking “install Hotfix 4.17.1.3-3.bat”
8. Follow instructions displayed in the window “Installation Cedex 2 Hotfix 4.17.1.3-3”

Subsequent to the installation of this hotfix the reagent kit has to be changed. This change has to be confirmed in the Cedex 2 software.

De-Installation

It is not possible to de-install the hotfix.

Validation

This hotfix does not influence the analysis process and has no impact on measured and calculated results!

Cedex 2 **HOTFIX**



Test

The installation of this hotfix was successfully tested with the following Cedex 2 software versions:

- Cedex 2.1.1 Build 4.16.1.0
- Cedex 2.1.1.Build 4.17.1.4
- Cedex 2.1.1 Build 4.20.1.0

Support

For further questions regarding the hotfix and its installation please contact:

support@innovatis.com

~ 082007.pm/ohe ~